



BioChip-D (Order No. 08503)

cellasys GmbH
Illerstrasse 14
87758 Kronburg / Germany

www.cellasys.com
info@cellasys.com

R&D:
Ohmstrasse 8
80802 München / Germany

Tel.: +49 89 2000110-74
Fax: +49 89 2000110-76

Content

Content	1
General	2
Description	2
Schematic (top side)	3
Schematic (bottom side)	4
Pin configuration	5
Technical data	7
pH (PH)	7
Dissolved oxygen (O2)	7
Impedance (IDES)	7
Temperature (TEMP)	7
Intended use	8
Intended misuse	8
Liability / Copyright	9

General

Please check delivery for transport damage when unpacking.

Description

Multiparametric BioChip for measurement (Impedance, pO₂, pH and temperature) of cellular vitality and changes in bioimpedance on glass substrate for optical access via microscope.

Caution


Handle with care.

- Glass tends to break due to mechanical stress.




In general **BioChip-D** may only be used in combination with cellasys IMOLA-IVD by **qualified personnel** of a research or healthcare institution. Read the **IMOLA-IVD manual** thoroughly and carefully follow the instructions and guidelines provided.

Schematic (top side)

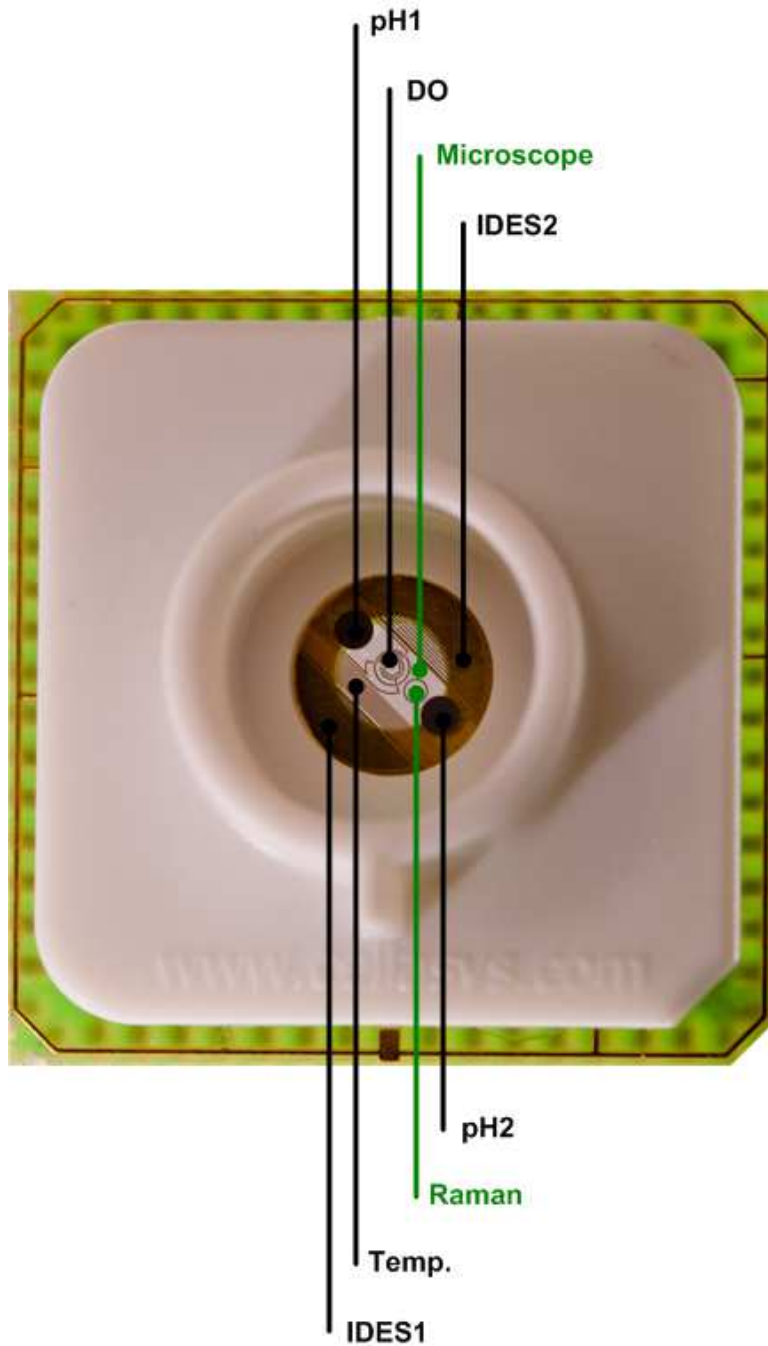

- IDES: Interdigitated electrode structure
- DO: Dissolved oxygen
- Temp: Temperature

Schematic (bottom side)

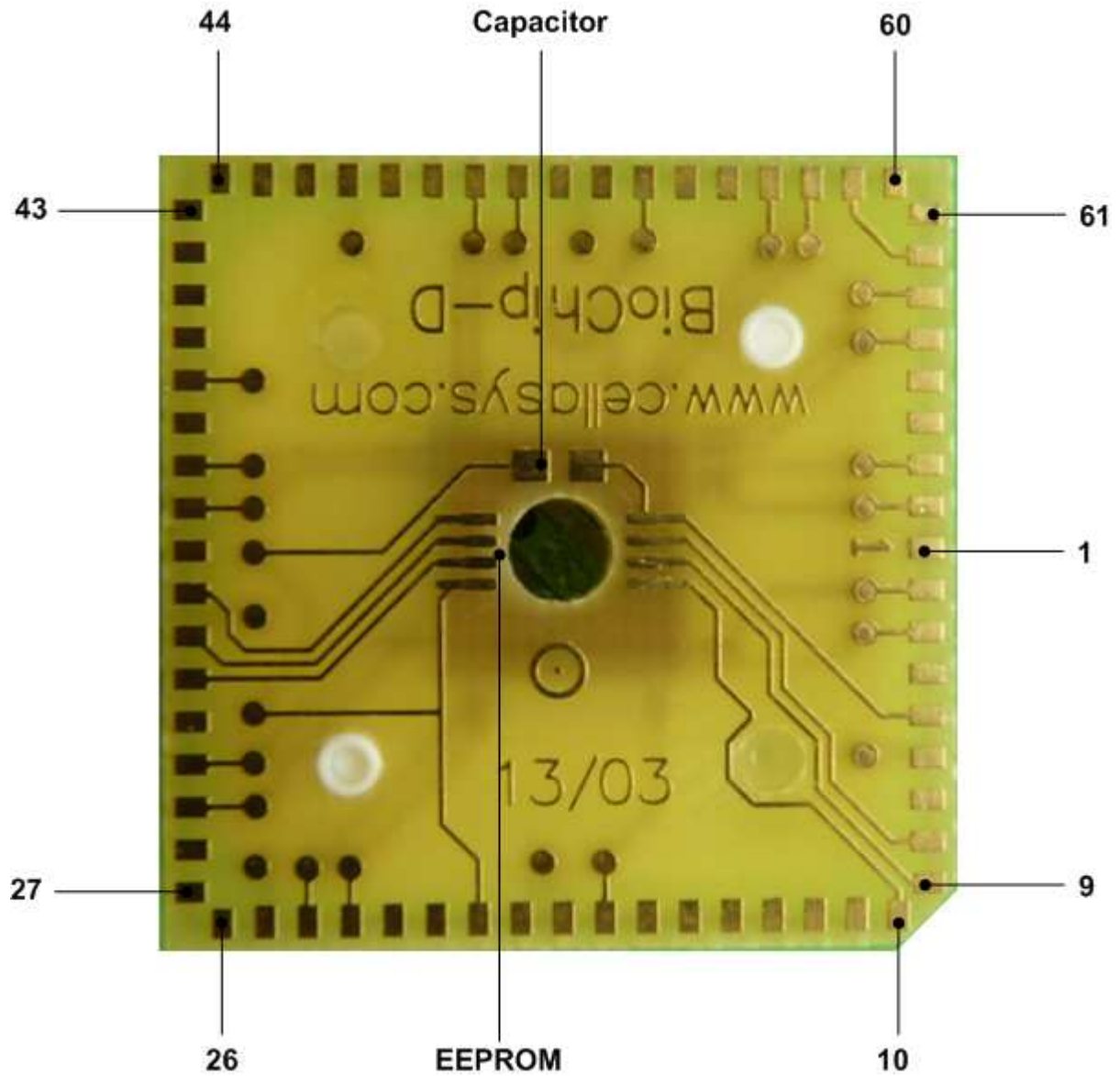

EEPROMs: M24C01 (TSSOP) – 401RP KTK726
 24AA08 (MSOP) – 4A08I 5123M1

Capacitor: 100nF (0603)

Pin configuration

Pin No.	Name	Description
1	-	Not connected
2	TEMPAU	Temperature sensor, voltage connector A
3	TEMPAI	Temperature sensor, current connector A
4	-	Not connected
5	MEMGND	EEPROM GND
6	-	Not connected
7	-	Not connected
8	MEME2	EEPROM chip enable 2
9	MEME1	EEPROM chip enable 1
10	MEME0	EEPROM chip enable 0
11	-	Not connected
12	-	Not connected
13	-	Not connected
14	-	Not connected
15	-	Not connected
16	-	Not connected
17	PH2	pH sensor 2
18	-	Not connected
19	-	Not connected
20	MEMVCC	EEPROM VCC (2.5V – 5.5V)
21	-	Not connected
22	-	Not connected
23	IDES2AU	Bioimpedance sensor 1, voltage connector A
24	IDES2AI	Bioimpedance sensor 1, current connector A
25	-	Not connected
26	-	Not connected
27	-	Not connected
28	-	Not connected
29	IDES2BI	Bioimpedance sensor 1, current connector B
30	IDES2BU	Bioimpedance sensor 1, voltage connector B
31	-	Not connected
32	MEMWC	EEPROM write control
33	MEMSCL	EEPROM serial clock
34	MEMSDA	EEPROM serial data
35	-	Not connected
36	O2REF	Dissolved oxygen sensor, reference electrode
37	PH1	pH sensor 1
38	-	Not connected
39	-	Not connected
40	-	Not connected

Pin No.	Name	Description
41	-	Not connected
42	-	Not connected
43	-	Not connected
44	-	Not connected
45	-	Not connected
46	-	Not connected
47	-	Not connected
48	-	Not connected
49	-	Not connected
50	O2WK	Dissolved oxygen sensor, work electrode
51	O2AUX	Dissolved oxygen sensor, auxiliary electrode
52	-	Not connected
53	-	Not connected
54	-	Not connected
55	-	Not connected
56	-	Not connected
57	IDES1AU	Bioimpedance sensor 2, voltage connector A
58	IDES1AI	Bioimpedance sensor 2, current connector A
59	SC	Shortened to 62
60	-	Not connected
61	-	Not connected
62	SC	Shortened to 59
63	IDES1BI	Bioimpedance sensor 2, current connector B
64	IDES1BU	Bioimpedance sensor 2, voltage connector B
65	-	Not connected
66	-	Not connected
67	TEMPBI	Temperature sensor, current connector B
68	TEMPBU	Temperature sensor, voltage connector B

Technical data

Dimensions: 24.0 x 24.0 x 10.0 mm³
 Weight: 4.5 g

Operating temperature: 0 °C to +80°C

pH (PH)

Dimensions (MeOx-Spot): ~ 3 mm²
 Linear range: pH 5.0 to pH 11.0
 Sensitivity: ~ - 40 mV/pH
 Response time (t₉₀): < 5 s

Dissolved oxygen (O2)

Dimensions: ~ 3 mm²
 Linear range: 0 to 120 %DO
 Sensitivity: 1 nA/pDO +/- 10 %
 Response time (t₉₀): < 0.1 s

Impedance (IDES)

Dimensions: ~ 10 mm²
 Linear range: 10 Ω to 5 kΩ
 Geometry: IDES1: 50 μm width, 25 μm distance
 IDES2: 50 μm width, 50 μm distance
 Response time (t₉₀): < 1 s

Temperature (TEMP)

Dimensions: ~ 4 mm²
 Linear range: 0 °C to +80 °C
 Sensitivity: 2.2 Ω/°C +/- 10 %
 Response time (t₉₀): < 1 s

Intended use

The BioChip-D is designed to be used in combination with IMOLA-IVD, for multiparametric measurement (impedance, pO₂, pH and temperature) of cellular vitality and bioimpedance.

The BioChip-D is a single-use device; it must not be used for multiple applications.

Intended misuse

The BioChip-D must not be operated with reagents and reagent products listed in directive 98/79 EEC, Annex II List A and List B.

Liability / Copyright

All technical details are state of the technology from February 2017 and are subject to change without notice. No liability is assumed for pictures, descriptions or any content of this document.

All descriptions, pictures, technical drawings and all other illustrations are protected by copyright and unless otherwise marked property of cellasys GmbH.

Any subsequent use needs prior written, allowance by cellasys GmbH.